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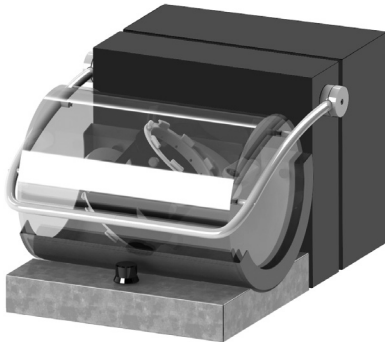
Data sheet Inversina 2 L (Tumbler mixer)

Bioengineering tumbler mixers are used to mix powders, solids, and liquids. Thanks to the efficient tumbling motion, constituents can be mixed quickly, gently, and thoroughly without the use of stirrers.

Bioengineering tumbler mixers are highly flexible as regards the deployment of mixing vessels. For instance, they can accommodate bottles, test tubes, or hoppers – pressurized, evacuated, heated, or cooled.

Interchangeable mixing vessels eliminate many cleaning processes as well as demixing caused by frequent decanting.

- Mixing of various solid and liquid substances.
- High mixing quality without the use of stirrers.
- No demixing of product as a result of frequent decanting, thanks to interchangeable mixing vessels.
- Deployment of interchangeable mixing vessels.
- Deployment of pressurized, evacuated, heated or cooled vessels.
- Easy to clean and maintain.

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Applications

The Inversina can be used to mix various types of substances.

Solid – solid

Solid-solid mixing is used mainly to mix powders for specimen formulation in analytical chemistry.

With appropriate abrasive media, the Inversina can also be used as a ball mill to debur components.

Solid – liquid

The solid-liquid combination is used to create slurries or to dissolve solids.

Liquid – liquid

Mixing liquids with liquids is a process used in paints, for example, for extraction purposes, or for producing emulsions.

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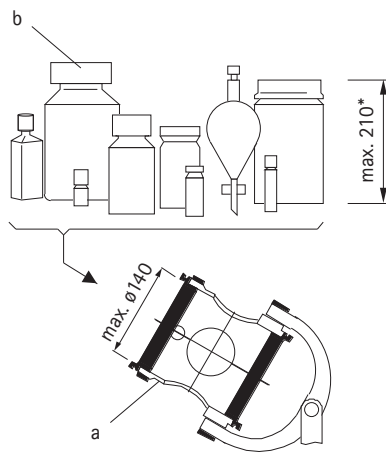
* The orientation of asymmetric vessels in the cage may reduce the maximum height; such vessels should be tested before being deployed.

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Product description

The Inversina is equipped with an electrically driven mixer cage (a) in which various mixing vessels (b) can be clamped.

The Inversina cage will accommodate mixing vessels (b) of various shapes with capacities of up to 2 L, maximum diameters of up to 140 mm, and maximum heights of about 210 mm. Mixing vessels such as wide-neck flasks can easily be secured in the cage by turning an rotating ring (a1) which tightens the clamping rings (a3).

The speed of the mixer cage can be adjusted at the control panel of the Inversina.

The transparent shroud (c) prevents inadvertent contact with moving parts. When it is open, the Inversina cannot be started; the Inversina also stops automatically, when the shroud is opened.

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